

The SC-FAST system for ICP and ICMS:

Versatile automation for routine and advanced environmental applications





FSEA Conference
May 28, 2010


Nathan Sævetit
Elemental Scientific




Elemental Scientific | esi@icpms.com | www.icpms.com

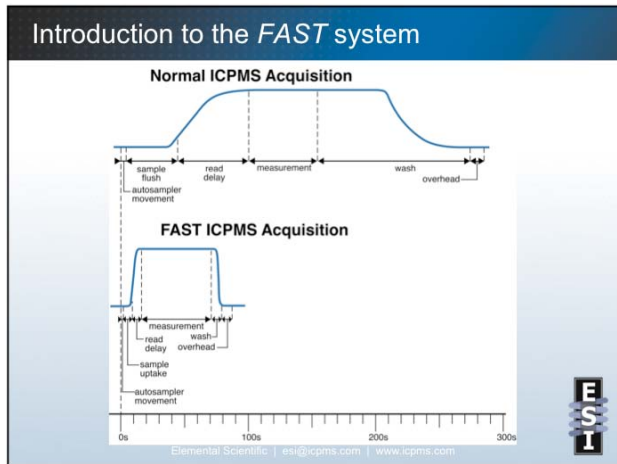

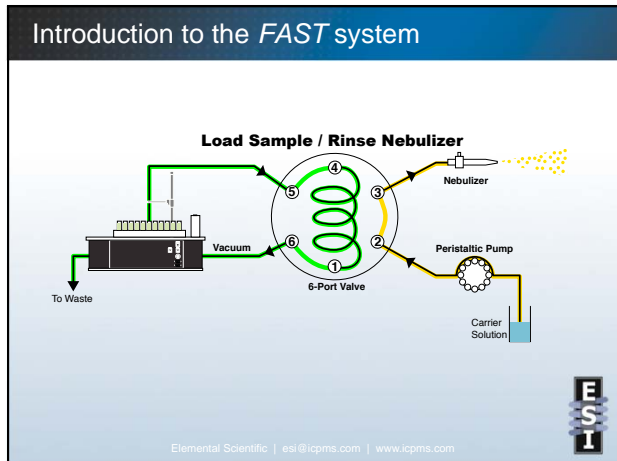

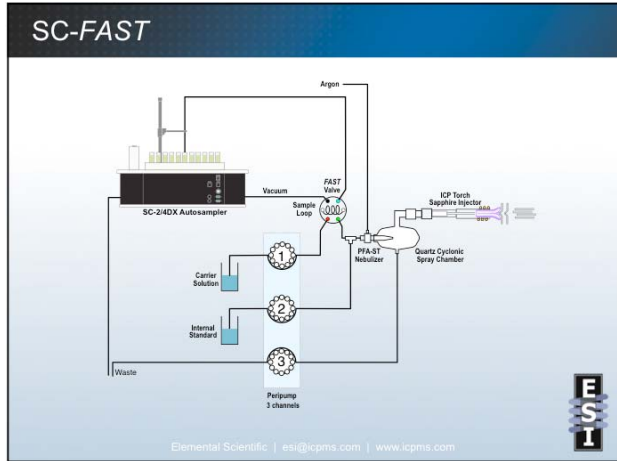

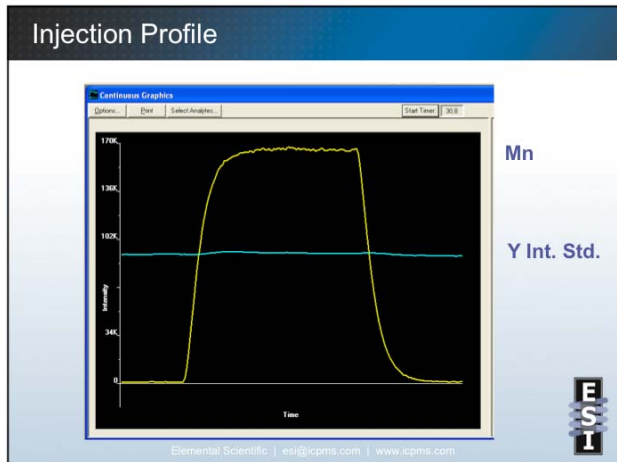

New SC-DX Autosamplers


- Enhanced chemical resistance
- Dual X-rails and heavy duty Z-rail
- Precise sampling, even on microtiter plates
- Optional enclosures and mobile stands improve cleanliness and convenience
- Easy integration of accessories

Elemental Scientific | esi@icpms.com | www.icpms.com

Selected FAST Applications

Advanced FAST Applications:


- seaFAST: Multimode analysis of undiluted seawater by ICPMS
- brineFAST: Multimode analysis of undiluted brines by ICPOES



Elemental Scientific | esi@icpms.com | www.icpms.com

seaFAST3


- seaFAST³ uses 3 modes of analysis for undiluted seawater or other high-matrix environmental waters:
 - Direct analysis with online dilution
 - Hydride generation
 - Preconcentration and matrix removal
- No sample preparation required
- Simple, external calibration with acid standards
- ~10 minutes, sample-to-sample
- All fluoropolymer sample path
- Full automation through ICPMS software




Elemental Scientific | esi@icpms.com | www.icpms.com

Modes for seaFAST3

H																	He		
Li	Be											B	C	N	O	F	Ne		
Na	Mg											Al	Si	P	S	Cl	Ar		
K	Ca	Sc	Ti	V	Cr	Mn	Fe	Co	Ni	Cu	Zn	Ga	Ge	As	Se	Br	Kr		
Rb	Sr	Y	Zr	Nb	Mo	Tc	Ru	Rh	Pd	Ag	Cd	In	Sn	Sb	Te	I	Xe		
Cs	Ba	Hf	Ta	W	Re	Os	Ir	Pt	Au	Hg	Tl	Pb	Bi	Po	At	Rn			
Fr	Ra																		
		La	Ce	Pr	Nd	Pm	Sm	Eu	Gd	Tb	Dy	Ho	Er	Tm	Yb	Lu			
		Ac	Th	Pa	U														






Elemental Scientific | esi@icpms.com | www.icpms.com


1) Direct Mode

- Online dilution
 - Reduces matrix effects
 - Reduces procedural blanks
- Online addition of internal standard
- Low-ppt calibrations




Elemental Scientific | esi@icpms.com | www.icpms.com

1) Direct Mode

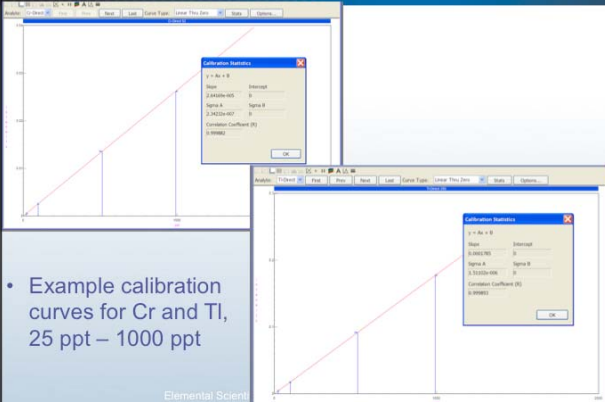


- Example calibration curves for Cr and Tl, 25 ppt – 1000 ppt

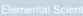


Elemental Scientific


2) Hydride Mode

- Specially designed batch hydride generation chamber
- High sample transport efficiency
- High sensitivity for As, Se, and Sb
- High purity syringe pumps automatically deliver the needed solutions at the required flow rates



Elemental Scientific | esi@icpms.com | www.icpms.com

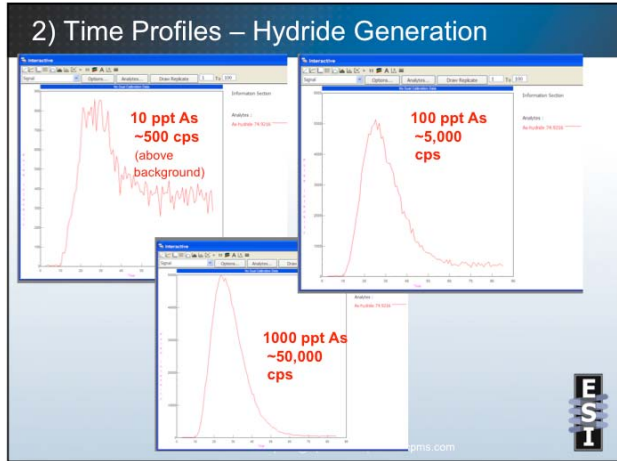

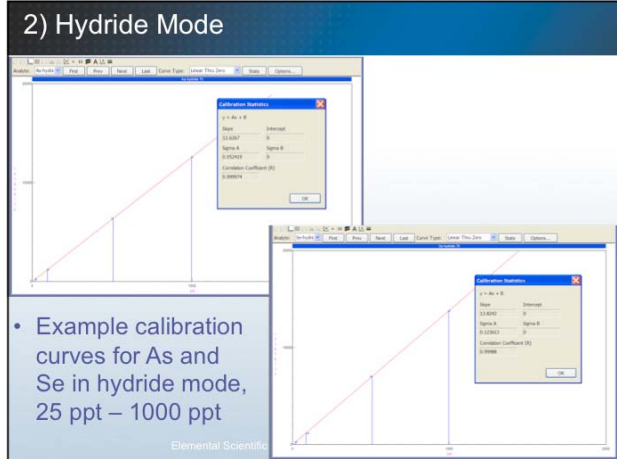



- Example calibration curves for As and Se in hydride mode, 25 ppt – 1000 ppt


3) Preconcentration and matrix removal

- Mixed resin containing iminodiacetate (IDA) resin
- Low affinity for alkali metal cations, alkali earth metal cations, and anions e.g. Na⁺, Mg²⁺, Ca²⁺, and Cl⁻
- High Affinity for many transition metal cations, M⁺

The image shows a white, cylindrical preconcentration column. It is labeled 'seFAST Concentrator' and 'High Purity PFA Column'. The ESI logo is in the bottom right corner.

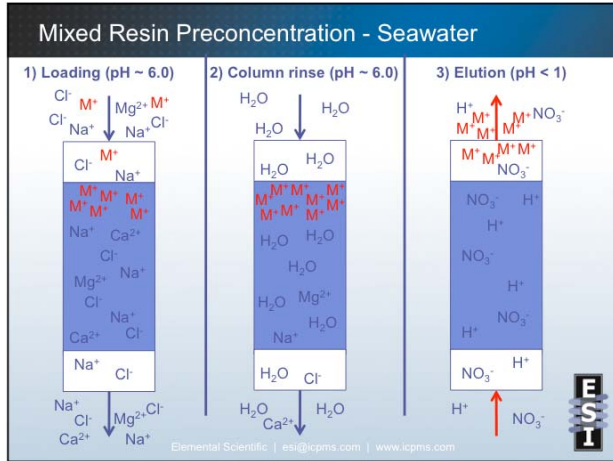

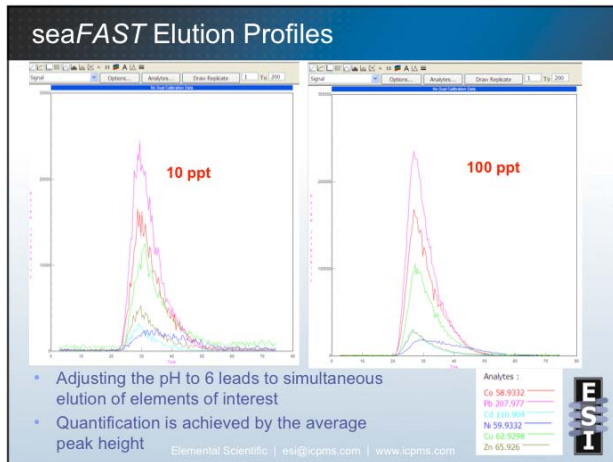

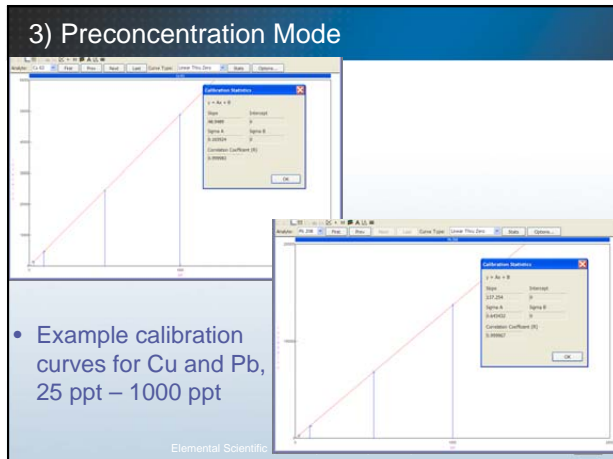


seaFAST3 - NASS-5 Recoveries with External Calibration

		Measured by seaFAST3	
		NASS-5	Certified (Reference)
Direct	Ag	312	n/a
	Ba	5382	(5100)
	Tl	12	n/a
Hydride	Se	14	(18)
	As	1146	1270 ± 120
	Sb	142	n/a
Preconcentration	V	1290	(1200)
	Mn	978	919 ± 57
	Fe	212	207 ± 35
	Co	12	11 ± 3
	Ni	259	253 ± 28
	Cu	287	297 ± 46
	Zn	183	102 ± 39
	Cd	25	23 ± 3
	Pb	7	8 ± 5


PerkinElmer Elan DRC
Elemental Scientific | esi@icpms.com | www.icpms.com




Method Detection Limits

Mode	Analyte	MDL, 25 ppt (ppt)
Direct	Ag	5
	Ba	8
	Tl	2
Hydride	Se	7
	As	3
	Sb	2
Preconcentration	V	2
	Mn	6
	Fe*	21
	Co	2
	Ni	7
	Cu	3
	Zn*	23
	Cd	3
	Pb	3

* Elements measured at 250 ppt, PerkinElmer Elan DRC
Elemental Scientific | esi@icpms.com | www.icpms.com



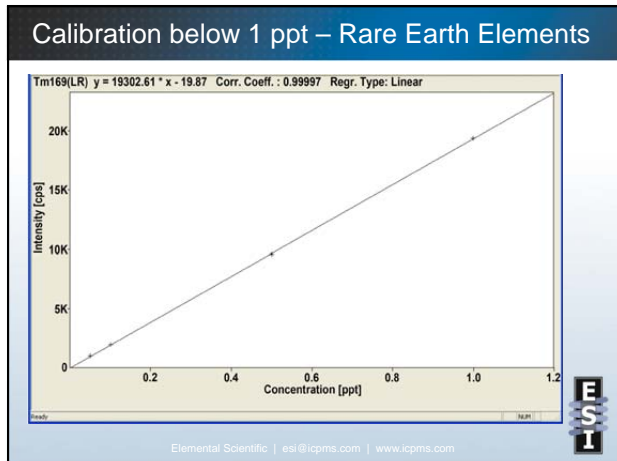


Rare Earth Element Determination with seaFAST

HRICPMS

Detection Limits (3σ) for Rare Earth Elements		Comparison of reference NASS-5 results and ESI seaFAST results (ppt)		
Analyte	LOD (ppq)	Analyte	Reference ¹	seaFAST
Y	12	Y	22.00 ± 2.90	24.17
La	29	La	12.19 ± 1.14	12.55
Ce	16	Ce	5.72 ± 0.69	5.93
Pr	5	Pr	2.08 ± 0.19	2.13
Nd	21	Nd	8.43 ± 0.65	8.80
Sm	11	Sm	4.74 ± 0.34	4.54
Eu	4	Eu	0.33 ± 0.03	0.32
Gd	6	Gd	1.83 ± 0.15	1.95
Tb	3	Tb	0.27 ± 0.03	0.25
Dy	6	Dy	1.62 ± 0.20	1.88
Ho	5	Ho	0.47 ± 0.06	0.42
Er	7	Er	1.43 ± 0.20	1.40
Tm	3	Tm	0.21 ± 0.04	0.20
Lu	4	Lu	1.29 ± 0.38	1.14
Yb	5	Yb	0.19 ± 0.06	0.20

Thermo ELEMENT2 HRICPMS
Elemental Scientific | esi@icpms.com | www.icpms.com


¹M. G. Lawrence and B. G. Saxe, *Geochimica et Cosmochimica Acta*, 2007, 71, 99-103.



brineFAST for ICPOES

- Similar chemistry to the seaFAST, except with a column that has retention for Ca and Mg, along with many other metals
- Ideal for ChlorAlkali plants or applications where Mg and Ca are analytes in high matrix samples
- Analyze undiluted, 30% brine without sample preparation
- Analyze caustic soda or bleach with minor sample preparation.


Elemental Scientific | esi@icpms.com | www.icpms.com



Calibration

- Simple calibration in acid standards
 - Low ppb calibration for ultrapure brines
 - May calibrate higher, 1ppm or more for more concentrated samples
- Same response in acid and brine
 - Spectral overlays
 - Spike recoveries

Elemental Scientific | esi@icpms.com | www.icpms.com




Method Detection Limits Comparison		
Analyte	Traditional (ppb)	brineFAST (ppb)
Mg	1.7	0.2
Ca	8.0	0.4
Mn	1.2	0.2
Fe	6.3	0.3
Sr	1.1	0.1
Ba	1.4	0.3
Al		2.0
Si*		8.6
Ti*		0.7
V		2.6
Cr*		2.1
Co		0.8
Ni		1.0
Cu		0.7
Zn		0.3
Cd		0.3
Pb		3.2



SC-FAST for ICP and ICPMS

- **Ideal for both high throughput labs and labs needing advanced automation or customization**
- Work on any ICP or ICPMS instrument
- Reduce operational costs
- Completely customizable
- Application solutions are available for advanced automation and flexibility
- Advanced online chemistries, hardware, and software overcome matrix effects
- Hardware and software can be configured for other applications as needed
- **Turnkey solutions for a wide variety of environmental analyses**

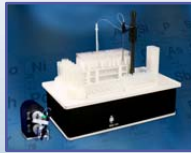

Elemental Scientific | esi@icpms.com | www.icpms.com
